



**A&A BIOTECHNOLOGY**  
innovating life science

## *Manual*

# Plasmid Maxi AX Sil

Increased efficiency kit for low- and high-copy plasmid DNA purification.  
Procedure with DNA precipitation.

catalog #	size
093-02S	2 isolations

For research use only.

### **Guarantee**

A&A Biotechnology provides a guarantee on this product.

The company does not guarantee the correct performance of this kit in the event of:

- not adhering to the supplied protocol
- use of not recommended equipment or materials
- use of other reagents than recommended or which are not a component of the product
- use of expired or improperly stored product or its components

# Table of Contents

<b>Specification</b>	<b>3</b>
<b>Contents</b>	<b>3</b>
<b>Additional equipment and reagents</b>	<b>3</b>
Necessary	3
Optional	3
<b>Important notes</b>	<b>4</b>
<b>Protocol</b>	<b>4</b>
<b>LySee color system</b>	<b>6</b>
Resuspension and lysis	6
Neutralization and precipitation	6
<b>Safety Information</b>	<b>7</b>

# Specification

<b>form</b>	maxicolumn
<b>binding capacity</b>	1 mg of DNA
<b>sample size</b>	up to 500 ml of bacteria culture
<b>elution volume</b>	precipitation

# Contents

<b>component</b>	<b>size</b>	<b>storage</b>
Maxi AX Sil 1000 columns	2 pcs	15–25 °C
Tubes 50 ml	6 pcs	15–25 °C
Filtration syringes	2 pcs	15–25 °C
L1 cell suspension solution	45 ml	2–8 °C
L2 lysis solution	45 ml	15–25 °C
L3 neutralizing solution	45 ml	15–25 °C
K1 Sil equilibrating solution	55 ml	15–25 °C
K2 Sil wash solution	130 ml	15–25 °C
K3 elution solution	55 ml	15–25 °C
TE buffer	5 ml	15–25 °C
Isopropanol	50 ml	15–25 °C

# Additional equipment and reagents

## Necessary

- Ethanol 70%
- Centrifuge
- Laboratory beakers

## Optional

- Sterile water (nuclease free) (cat.# 003-075, 003-25)

## Important notes

- Kit contains the LySee color system for easy optical control of alkaline lysis progress (page 6).
- SDS detergent is a component of L2 lysis solution and precipitates at low temperatures. Whenever the L2 lysis solution is not clearly transparent it must be warmed at 40 °C to form a thoroughly clear solution.

## Protocol

1. Centrifuge **200 to 500 ml** of overnight bacterial culture.

**Note.** Volume of bacterial culture that can be processed depends on the particular plasmid copy number. Please keep in mind that the total binding capacity of the column is 1 mg DNA.

2. Discard supernatants. Suspend the bacterial pellet in **20 ml** of **L1** cell suspension solution.

**Note.** During the bacterial pellet suspension, the solution will change color from a transparent deep pink to opaque light pink. Suspending is considered completed when the pellet, at the bottom of the tube disappears.

3. Prepare two **tubes 50 ml** (included). Transfer **10 ml** of the suspension to each of the two tubes.

4. Add **10 ml** of **L2** lysis solution to each tube. Gently mix by inverting the tubes.

**Note.** After the addition of L2 lysis solution, gently mix the tube contents so as not to cause fragmentation of the chromosomal DNA. Mix the tube by inverting (5-6 times). The mixture should change appearance and color.

5. Keep for **3 min** at **room temp.**

**Note.** After 3 min, the lysate should be clear and raspberry in color. If not, mix the lysate several times and leave for an additional 3 min.

6. Add **10 ml** of **L3** neutralizing solution to each tube. Gently mix until the raspberry color disappears.

**Uwaga.** After the addition of L3 neutralizing buffer followed by the rapid precipitation of the potassium salts (SDS), chromosomal DNA and certain proteins. After mixing, the tube contents should change the color to yellowish. No traces of raspberry color indicates complete neutralization and successful ending of the alkaline lysis.

7. Keep on ice for **15 min.**

8. Centrifuge for **10 min** at **10 000 RPM.**

9. Prepare the Maxi AX Sil 1000 column by placing it in the upper opening of the rack. Place a beaker under the column to collect waste.

Apply **25 ml** of **K1 Sil** equilibrating solution onto the Maxi AX Sil 1000 column. Wait for the solution to flow through the column.

10. Remove the plunger from the **filtration syringe** and pour the supernatants (with leftovers of precipitate) of both 50 ml tubes (point 8.) into the **filtration syringe.**

11. Insert the plunger and filter the lysate directly to the Maxi AX Sil 1000 column. Wait for the lysate to flow through the column.
12. Add **30 ml** of **K2 Sil** wash solution to the Maxi AX Sil 1000 column. Wait for the solution to flow through the column.
13. Again, add **30 ml** of **K2 Sil** wash solution to the Maxi AX Sil 1000 column. Wait for the solution to flow through the column.
14. Remove the beaker under the column Maxi AX Sil 1000. Place a **new 50 ml** tube (included) under the column.
15. Add **25 ml** of **K3** elution solution to the Maxi AX Sil 1000 column. Wait for the eluate to flow through the column.
16. Remove the column Maxi AX Sil 1000 from the rack. Add **20 ml** of **isopropanol** to the tube.
17. Close the tube, mix and centrifuge for **20 min** at **11 000 x g**.
18. Carefully discard supernatant. Be careful not to remove the DNA pellet at the bottom of the tube.  
**Attention!** When pouring out the supernatant, it is very easy to lose the DNA pellet. To be safe, it is recommended to pour the supernatant into the prepared tube so the pellet can be recovered.
19. Add **2 ml** of **70% ethanol** (not included). Gently mix and centrifuge for **3 min** at **11 000 x g**.
20. Carefully discard supernatant. Be careful not to remove the DNA pellet at the bottom of the tube.  
**Attention!** When pouring out the supernatant, it is very easy to lose the DNA pellet. To be safe, it is recommended to pour the supernatant into the prepared tube so the pellet can be recovered.
21. Air dry the plasmid DNA pellet for **10 min** at **room temp.** up-site down.  
**Note.** If there are any leftovers (small droplets) of alcohol on the tube walls they should be removed with sterile cotton buds.
22. Dried DNA pellets can be dissolved in **0.2-1 ml** of **TE** buffer or **sterile water** (not included). Store the plasmid DNA at 4-8 °C.

## LySee color system

The LySee color system enables an easy and convenient visual control of alkaline lysis. The visual control system prevents common handling errors of incomplete cell resuspension, inefficient cell lysis and incomplete precipitation of unwanted cell components.

### Resuspension and lysis

The addition of the transparent purple L1 color cell suspension solution to the bacterial cell pellet makes the bacterial cell pellet easy to localize (fig 1). During the suspension of the bacterial cell pellet, the solution turns opaque light pink (fig 2). The suspension is completed with the complete disappearance of the pellet at the bottom of the tube. After the addition of L2 lysis solution and incubation, lysate turns transparent raspberry. Cell lysis is completed when the solution will turn homogeneously transparent raspberry (fig 3).

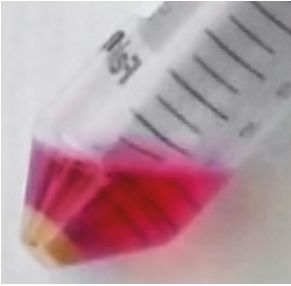


fig 1

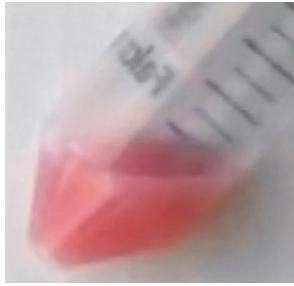


fig 2

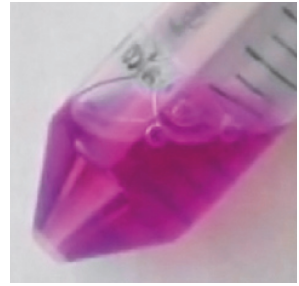


fig 3

### Neutralization and precipitation

The addition of the L3 neutralizing solution causes rapid precipitation of potassium salts (SDS), chromosomal DNA and some proteins (fig 4). After mixing, the solution turns yellowish (fig 5). No traces of raspberry color indicates complete neutralization and successful ending of alkaline lysis (fig 6).

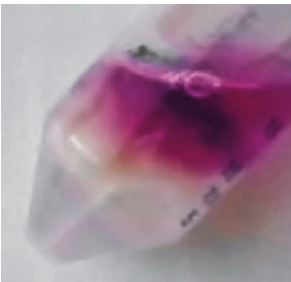


fig 4



fig 5



fig 6

# Safety Information

---



**WARNING**

## L2 lysis solution

H315 Causes skin irritation.  
 H319 Causes serious eye irritation.  
 H334 May cause allergy or asthma symptoms or breathing difficulties if inhaled.  
 H335 May cause respiratory irritation.  
 P261 Avoid breathing dust.  
 P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.  
 P342+P311 If experiencing respiratory symptoms call a Poison Center or doctor/physician.

---



**DANGER**

## K3 elution solution

H225 Highly flammable liquid and vapor.  
 H319 Causes serious eye irritation.  
 H336 May cause drowsiness or dizziness.  
 P210 Keep away from heat, sparks, open flames, hot surfaces. No smoking.  
 P261 Avoid breathing vapors.  
 P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

---



**DANGER**

## K1 Sil equilibrating solution

H225 Highly flammable liquid and vapor.  
 H319 Causes serious eye irritation.  
 H336 May cause drowsiness or dizziness.  
 P210 Keep away from heat, sparks, open flames, hot surfaces. No smoking.  
 P261 Avoid breathing vapors.  
 P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

---



**DANGER**

## K2 Sil wash solution

H225 Highly flammable liquid and vapor.  
 H319 Causes serious eye irritation.  
 H336 May cause drowsiness or dizziness.  
 P210 Keep away from heat, sparks, open flames, hot surfaces. No smoking.  
 P261 Avoid breathing vapors.  
 P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

---



**DANGER**

## Isopropanol

H225 Highly flammable liquid and vapor.  
 H319 Causes serious eye irritation.  
 H336 May cause drowsiness or dizziness.  
 P210 Keep away from heat, sparks, open flames, hot surfaces. No smoking.  
 P261 Avoid breathing vapors.  
 P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

---



**A&A BIOTECHNOLOGY**  
innovating life science

A&A Biotechnology, ul. Strzelca 40, 80-299 Gdańsk, Poland  
phone +48 883 323 761, +48 600 776 268  
info@aabiotech.com, www.aabiotech.com

